**AUTOCLAVI  
760-770**

**AUTOCLAVES  
760-770**

MADE IN ITALY

Distribuito da:  
Zetalab s.r.l.  
Via Castelfidardo, 11 - 35141 Padova  
Telefono 049 2021144 - Fax 049 2021143  
Internet: [www.zetalab.it](http://www.zetalab.it) - e-mail: [info@zetalab.it](mailto:info@zetalab.it)

**Zetalab**

---

## AUTOCLAVE “VAPOR MATIC” 770

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---

## AUTOCLAVE “VAPOR MATIC” 770

The Vapor Matic 770 is a sterilization autoclave with vertical charging which is completely automatic, thermoregulated and microprocessor controlled. The Vapor Matic 770 is equipped with a microprocessor with high disturbance resistance HMOS logic, which permits the programming of various times and temperatures, including the two most common sterilization cycles: cycle 1 = 121 °C – 1 Atm – 20'; cycle 2 = 134 °C 2 Atm – 20'. These cycles can be memorized and called up by the pressing of a single key so the setting of the autoclave is immediate. F.U. cycle switch 30-60'.

After the starting of the process a valve eliminates the air, present in the boiler and the initial condensation. After the programmed temperature is reached this valve automatically closes and the display gives the count-down of the time required to complete the process. The end of the process is signaled by an audible signal.

**In case of temporary power failure, the sterilization is automatically restored exactly from the point it was interrupted. In long power cut-offs of there is a 3 °C difference from preset temperature rate, the full sterilization time is automatically set and the whole cycle is carried out again.**

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### Technical specifications

- Outer casing in powdered painted.
- Inner camera in aisi 304.
- Shielded heating element located al the bottom.
- Protection lack of water.
- Cover with 1 knobs and silicone gasket and protection cover.
- Cover security system.
- Signaling lamp when temperature >80 °C.
- Operating temperature: 100 °C-134 °C (max. pressure 2.0 Atm).
- Temperature registering probes PT 100.
- Safety valves set 2.5 Atm.
- Air and condensed relief valve.
- Cock for complete water drain from the boiler.
- Manometer for pressure indication.
- Sterilization time selectable from 1 up to 99 minutes.
- Power supply: 230 V – 50 Hz – 2200 W.
- Weight: Kg.63
- Not subject to I.S.P.E.S.L. testing being within capacity and pressure limits.

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### Error codes

- *Safety self-diagnosis which advises the operator by means of audible and visible warnings (danger led) of an incorrect temperature setting (=142 °C) or an internal fault in the power circuit.*
- *Alarm signalling at the temperature Display Unit for:*
  - lid open;
  - no water;
  - PT 100 circuit open;
  - PT 100 short circuit.
  - Temporary power failure.

### Accessories

- 2 closed stainless steel baskets (Ø 240 x 190 mm. ht.).
- 1 perforated stainless steel basket (Ø 240 x 430 mm. ht.)
- 500 pieces packet of heat proof bags (360 x 750 mm. ht.).
- Electronic printer extra strong paper.

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
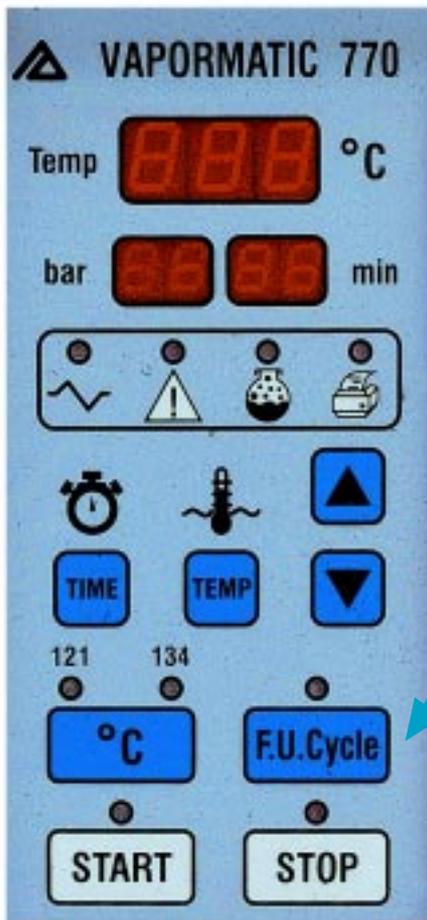
### OFFICIAL PHARMACOPEIA CYCLE to sterilize as follow:

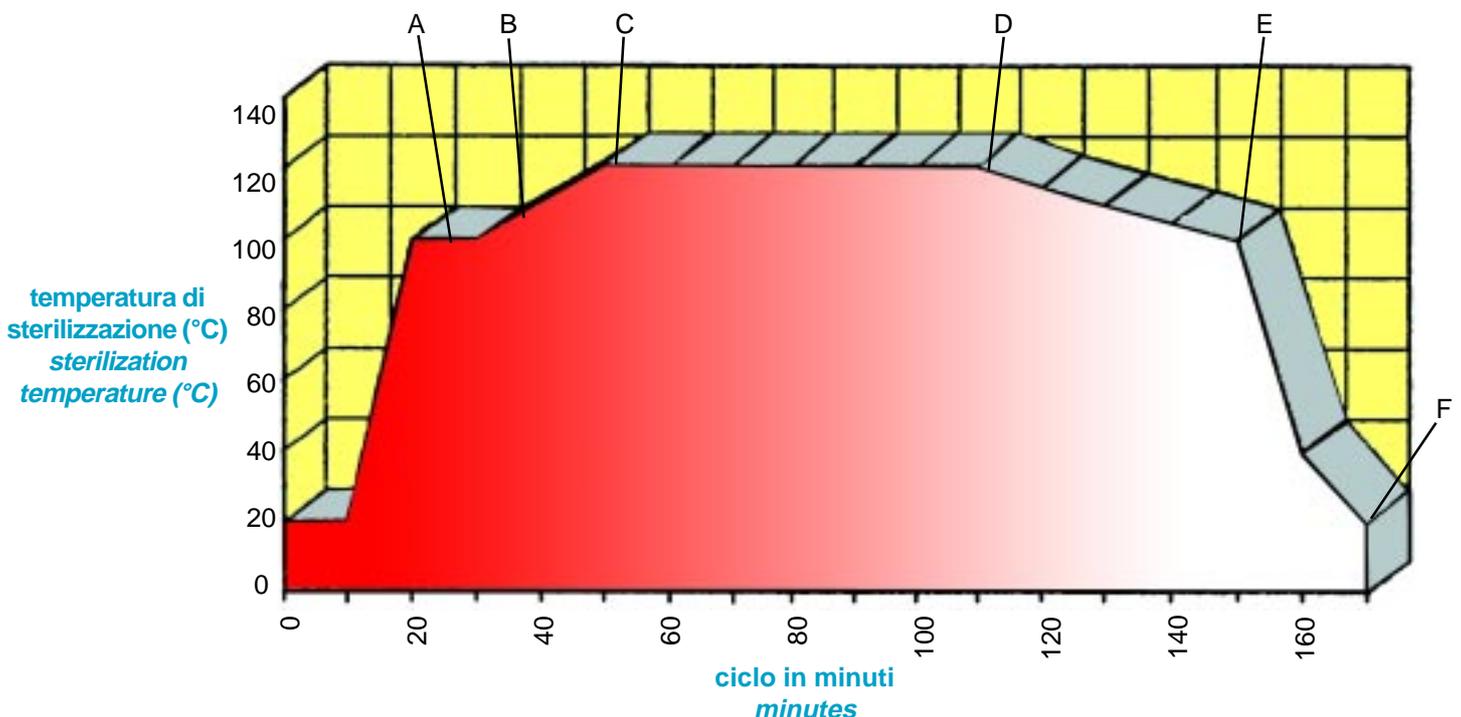- Up to 100°C and stabilization for 10' to eliminate air (A)
- Up to 121°C for 1°C/min. total 20' (B)
- Sterilization for 30'/60' at 121°C (1 Atm) (C)
- Down temperature at speed 0.5°C/min. (D)
- End down at 100°C in 40'.(E)
- Down temperature until room temperature (F)

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## AUTOCLAVE 760

*Sterilization chamber in aluminium. Lid with fastener fitted with safety and exhaust valves.*

*Manometer for working pressure indication; radiant electric heating plate.*

*Protection device for insufficient level of the water in the tank catches fire the spy "lack water".*

*Maximum working pressure: 1 Atm (121 °C). Outer finish in powdered painted steel*

- Power supply: 230 V – 50 Hz – 1500 W.
- Weight: 18 Kg.

*Not subject to I.S.P.E.S.L. testing being within capacity and pressure limits.*

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## Accessories

- 1 closed stainless steel basket (Ø mm. 240 x 240ht.)
- 1 perforated stainless steel basket (Ø mm. 240 x 270ht.)
- 500 pieces packet of heat proof bags (mm. 360 x 750ht.)

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**CE** Is in conformity with the following Directives: CEE – 73/23, 89/392, 89/336, 91/368, 92/31, 93/44, 93/68 the following Rules: CEI EN 61010- 1 (IEC 1010-1/CEI 66-5) - the following National Laws: D.P.R. N. 547 27/4/1955 Law N. 186 1/3/1968