## GC ZERO AIR GENERATOR

The GC Zero Air generators produce dry and hydrocarbon-free air, using air from oil-free compressors, thus avoiding the need to use traditional bottles that are often complex to change.

The ergonomic, compact and silent design makes the GC series generators the ideal solution for the following analytical applications: GC-FID, GC-NPD, GC-FPD, GC-PFPD, THA.

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#### EXTERNAL COMPRESSOR (option)

Following an agreement with a leading manufacturer operating a sales and service network in many cities throughout Italy and the world, DBS can supply the right "oil free" compressor for different flow requirements.

### OPERATING DIAGRAM

Zero Air generators use three steps to transform ambient air into analytical grade air.

#### STEP 1: PRE-FILTRATION

The external oil-free compressor delivers air through a high efficiency filter that traps any particles that may damage the system. The filter has an automatic purge system, and traps oil, water and any other particles larger than 5 microns in size.

#### STEP 2: HC AND CO TRAPPING

The air leaving the filter enters a high-temperature platinum catalyser, which through oxidation eliminates hydrocarbon molecules down to < 0.1 ppm.

#### STEP 3: FINAL FILTRATION

A high efficiency filter is used to prevent any kind of particles from entering the instrument.

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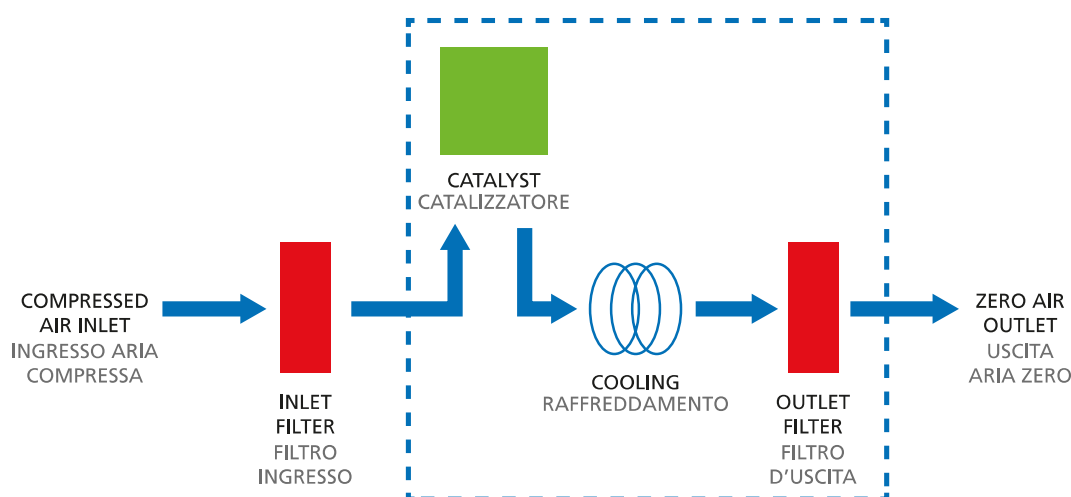
**BENEFITS****BETTER DETECTOR PERFORMANCE**

The reduction of hydrocarbons, including methane, and carbon monoxide to < 0.1 ppm decreases the background noise level and gives the baseline much better stability, considerably increasing detector sensitivity and ensuring precise analytical results.

**EASY INSTALLATION**

Just connect a suitable oil free compressor (optional) to almost immediately have air at the right grade.

To save bench space, the unit can be easily installed on the lab wall.

**LAB SAFETY**

No more bottles in the lab and expensive pipelines for air distribution.

**SAVE MONEY**

The unit only requires connection to a suitable oil free compressor and the mains: the investment can be paid back in less than a year.

**IMPROVED LAB EFFICIENCY**

Continuous operation 24 hours a day cuts dead times for gas bottle changeover and avoids the need for tedious instrument recalibrations.

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