MINI WHISPER

The Mini-Whisper nitrogen generator has been developed to meet specific requirements in terms of flow, purity and pressure on some LCMS interfaces as well as on light scattering evaporative detector.

The simple high efficiency membrane technology allows the separation of nitrogen from the other components of the compressed air inlet.

The low pressure drop allows the unit to be connected to an existing dry and oil-free compressed air source in the lab.

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None

ROUTINE MAINTENANCE

Simple filter replacement once a year.

BENEFITS

BETTER RESULTS FROM ANALYTICAL INSTRUMENTS

The constant purity of the nitrogen improves system stability and ensures reproducible results.

BETTER LAB EFFICIENCY

Gas bottles are insufficient to ensure good lab productivity.

COST SAVINGS

The investment can normally be paid back in less than a year. No pipeline is required to carry the gas from the storeroom to the lab.

IMPROVED SAFETY

Nitrogen produced at low pressure and at room temperature eliminates the risks relating to high pressure gas bottle use and the handling of liquid nitrogen.

...AND SIMPLE INSTALLATION

The Mini-Whisper nitrogen generator is very easy to install in the lab. No power connection is required. Bench-top or wall-mounted installation to save bench space!

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**MEMBRANE SEPARATION TECHNIQUE**

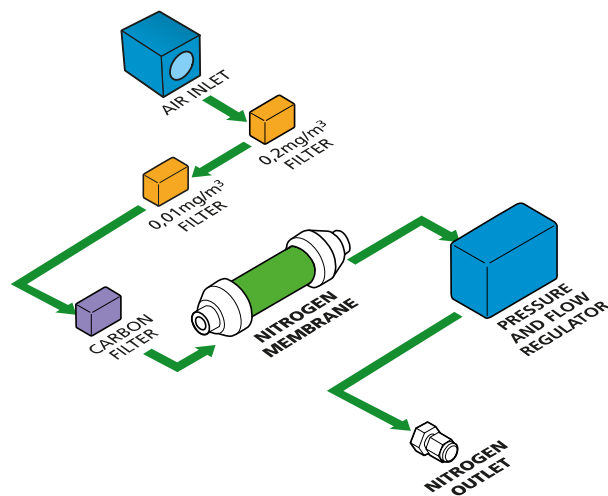
Compressed air is forced through a hollow fibre membrane, with selective permeation of the different components - nitrogen, oxygen, CO₂, water vapour and traces of rare gases - depending on the rate of diffusion.

The rate of diffusion of nitrogen through the membrane is slower than the other components, which thus flow outside of the fibre membrane, leaving just nitrogen inside; this is then discharged subsequently, ready for use.

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MODELS modelli	N2-MINI-WHISPER
Out max flow	8-12 NI/min
Input air flow	> 40 lit/min
Out pressure	7 bar
Input air pressure	5-8 bar
Pressure drop	< 0.8 bar (10 psig)
Purity (based on outlet flow)	> 98 - 99.5%
Residual particulate	< 0.01 micron
Dewpoint	-40°C
Maximum operating temp	10°C – 35°C
Air compressor supplied	NO (optional)
Background noise	none
Electrical requirements	none
Connections	¼ G
Weight (kg)	8
Dimensions (mm)	735H x 348W x 350D