BILANCIA IDROSTATICA ELETTRONICA ELECTRONIC HYDROSTATIC BALANCE

SUPERALCOMAT

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ELECTRONIC HYDROSTATIC BALANCE
Fully automatic to determine Alcoholic
Strenght by Volume ASV, total dry
extract and musts evaluation.
Commission Regulation (EC) 2870/2000
and 128/2004.

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The readout of the Alcoholic Strenght by Volume is shown on a large LCD screen simultaneously with the readout of temperature, density at 20°C and other parameters.

The reading module is provided with a shower proof keyboards whose functional keys allow:

- correction of the floater volume with H₂O or certified standard solutions
- zeroset and calibration of the SUPER ALCOMAT
- date, hour, minute setting and other functions
- the printing of the reading values and other parameters (printer not included)
- musts evaluation: Baumé, Babo, Brix and Oechsle degrees
- working temperature range: 10 - 30°C optimum 15 - 25°C)
- languages: Italian, German, English, French and Spanish.

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A Motorola micro-controller, specially designed for Gibertini, besides managing Super Alcomat most important functions, it is able to detect any internal problem inside the balance that could affect its accuracy, and suggest a service solution carried out by our qualified staff.

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Readability

0,01 %ASV

Readability and reading repeatability

± 0,03 %ASV (between 15 and 25 °C)

TOTAL DRY EXTRACT

Range of values

0,0 ÷ 505,8 g/l of sucrose

Readability and reading repeatability

± 0,3 g/l

DENSITY

Measurement range for the relative density

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Standard equipment

Power pack

2 hydrorepellent floaters, interchangeable

in mass and volume

Container for exhaust liquids

Certified hydroalcoholic solution

Response time

6 sec.

Autocalibration with internal mass

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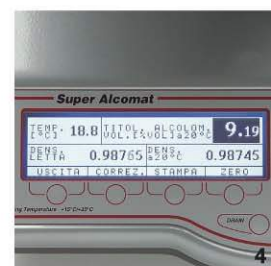
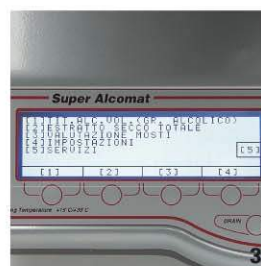
1. When the "drain" is open, flashing lamp is on 2. Automatic device to lift up and lower the floater 3. Main menu 4. Asv and density reading example 5. Device to lock the internal balance 6. Rear side: RS232 and printer outputs

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Floaters with certificate - SIT certified
solutions 5 ÷ 60 %ASV

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FLOATERS INTERCHANGEABLE IN MASS AND VOLUME

A particular care is devoted to the construction and calibration of the floaters.

They are all interchangeable in mass and volume to simplify their use; for example, it is possible to use two different floaters to determine the relative density of a must or a wine; in this way possible sample pollution is avoided.

In order to achieve what mentioned before, the suspension wire realised with a non-magnetic and stainless material, with a diameter less or equal to 0.20 mm. and the superficial treatment of the floater's glass were object of thorough studies to nullify the problems caused by the surface tension, by the wire's own volume, and by the electrostatic charges.

The floaters volume can, on request, be certified by competent organisations, such as the L.N.E. Paris, the I.M.G.C. Turin, the PTB Germany, etc.

Soluzioni idroalcoliche certificate SIT dal 5% al 60% v/v - a richiesta SIT Certified hydroalcoholic solutions from 5% to 60% v/v - on request

laboratory electronic balances & oenological equipments